Problem: Determine the density of water, and identify 4 metals based on their density.

Procedure:

WATER:

- 1. Find the mass of a clean, dry beaker with the digital balance (to the nearest cg or 0.01 g).
- 2. Fill the buret up to 0.00 (not past) with DISTILLED WATER and record the initial volume (nearest 0.01 mL).
- 3. Add up to 50 mL (not more) water from the buret into the empty beaker and record the final volume.
- 4. Find the mass of the beaker with water.

Mass clean dry beaker _____ g
Buret initial volume ____ mL
Buret final volume ____ mL

METALS:

- 5. Measure the mass of each metal cylinder with the digital balance (to the nearest cg or 0.01 g).
- 6. Find the diameter and height of each cylinder with the vernier caliper.
- 7. For one metal cylinder (since they are all similar) find the volume by water displacement (nearest mL). Take a PLASTIC graduated cylinder with enough water to submerge the metal. Read the initial water volume, add the metal and read the final water volume to find the metal's volume.
- 8. For the rectangular prism metal object, find the mass and dimensions (l,w,h) using the caliper.

Data:

Mass beaker + water g					
METAL					
Metal	#1	#2	#3	#4	prism
Color					
Mass (g)					
Diameter (mm)					L=mm
					W= mm
Height (mm)					
Water displacement:					
Initial volume	mL				
Final volume	mL				